

## HOW TO MOUNT POINT COUNTER (CAT. NO. MA945)

1. Attach the point counter to the rotating stage by inserting guide pins ① into the holes ② on rotating stage.
2. Tighten the set screw ③ by supplied Allen wrench ④. In case the set screw is hidden, move the point counter back or forward by turning counting knob ⑤ so that the set screw ③ coincides with female threads on the rotating stage.
3. When putting or changing counting knob ⑤, be sure to put it in order to engage the detent inside the knob cap.

NOTE: Figure on counting knob means travel length of the counter by one click movement.

4. Then, put specimen on the stage and hold it with finger ⑥.

